

# StepOnePlus™ Compatible Consumables


- 樣品量多時

- **P/N 4346907**

MicroAmp® Fast 96-Well Reaction Plate (0.1 mL) - 10 plates

- **P/N 4360954**

MicroAmp® Optical Adhesive Film - 25 films

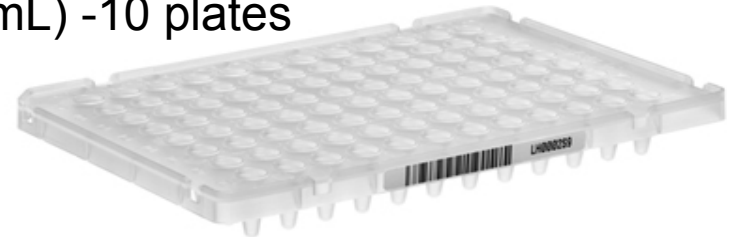


- 樣品量少時

- **P/N 4358293**

MicroAmp® Fast 8-Tube Strip (0.1 mL) - 125 strips

- **P/N 4323032**

MicroAmp® Optical 8-Cap Strip - 300 strips

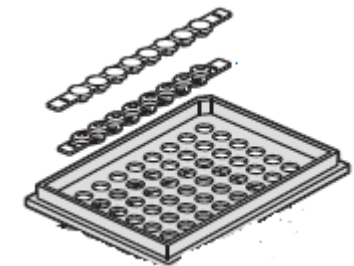

★ Place the tray containing the tube, Load at least 16 tube

## Operation Notes

- **Directly load fast optical 96-well plate into the instrument**
  - ✓ If using fast 8-tube stripes, load the tubes with fast 96-well tray
- Do not mark any labels on the consumables
  - ✓ This may increase the background signal
- Avoid bubbles when pipetting into each well
  - ✓ Centrifuge samples

